



**UV-4600**

**UV-VIS SPECTROPHOTOMETER**

# **USER'S MANUAL**

Version: August 2016

Mark



**DINTER**<sup>sa</sup>

Encarnació, 123 -125. Phone. +34 93 284 69 62. Fax +34 93 210 43 07

E-mail: [dinter@dinko.es](mailto:dinter@dinko.es) [www.dinko.es](http://www.dinko.es) 08024-BARCELONA

# Contents

<b>Chapter 1 Preface</b> .....	<b>4</b>
1. Safety.....	4
2. General.....	4
3. Electrical .....	4
4. Warning.....	4
5. Performance.....	5
6. Radio Interference.....	5
<b>Chapter 2 Introduction of the Instrument</b> .....	<b>6</b>
1. Brief Introduction of the instrument.....	6
2. Working Principle.....	7
3. Structure.....	7
4. Unpacking Instructions.....	8
5. Specifications.....	8
<b>Chapter 3 Install Instrument</b> .....	<b>10</b>
1. Environment Required.....	10
2. Voltage Confirmation.....	10
3. Installation.....	10
<b>Chapter 4 Operation Introduction</b> .....	<b>12</b>
1. Software Structure.....	12
2. Basic Operation.....	13
2.1 Choose Photometric Mode.....	13
2.2 Set Wavelength .....	13
2.3 Set Parameters.....	13
2.4 Set the position of auto-cell holder.....	13
2.5 Delete the Input Value.....	13
2.6 Delete the test result and stored data.....	13
2.7 Calibrate 100%T/0Abs.....	13
2.8 Measure Samples.....	13
2.9 Print the test result.....	14
3. Preparation before test.....	14
3.1 Switch on the instrument and begin the system self-test.....	14
3.2 Pre-warm.....	14
3.3 Check the cuvettes.....	14
4. Measurement.....	14
4.1 Basic Mode .....	16
4.2 Quantitative.....	16
4.3 Setup .....	23

1) Power On/Off D2 Lamp.....	23
2) Power On/Off W Lamp.....	23
3) Set Date and time.....	24
4) Calibrate Dark Current.....	24
5) Wavelength Calibration.....	25
6) Input the lamps changing Wavelength Point.....	25
7) Reload Default.....	26
8) Version.....	26
<b>Chapter 5 Instrument Maintenance.....</b>	<b>27</b>
1. Daily Maintenance.....	27
2. Trouble shooting.....	27
3. Spare part Replacement.....	28
<b>Accessories.....</b>	<b>29</b>
<b>CE Statement.....</b>	<b>30</b>

# CHAPTER 1 PREFACE

## 1, Safety

The safety statements in this manual comply with the requirements of the HEALTH AND SAFETY AT WORK ACT, 1974.

Read the following before installing and using the instrument and its accessories. This instrument should be operated by appropriate laboratory technicians.

## 2, General

The apparatus described in this manual is designed to be used by properly trained personnel in a suitable equipped laboratory. For the correct and safe use of this apparatus it is essential that laboratory personnel follow generally accepted safe procedures in addition to the safety precautions called for in this manual.

The covers on this instrument may be removed for servicing. However, the inside of the power supply unit is a hazardous area and its cover should not be removed under any circumstances. There are no serviceable components inside this power supply unit. For this instrument, avoid touching the high voltage power supply at all times.

Some of the chemicals used in spectrophotometry are corrosive and/or inflammable and samples may be radioactive, toxic, or potentially infective. Care should be taken to follow the normal laboratory procedures for handling chemicals and samples.

## 3, Electrical

Before switching on the apparatus, make sure it is set to the voltage of the local power supply (see **Installation**).

The power cord shall be inserted in a socket provided with a protective earth contact. The protective action must not be negated by the use of an extension cord without a protective conductor.

## 4, Warning

Any interruption of the protective conductor inside or outside the apparatus or disconnection of the protective earth terminal is likely to make the apparatus dangerous. Intentional interruption is prohibited.

Whenever it is likely that the protection has been impaired, the apparatus shall be made inoperative and be secured against any unintended operation.

**Note: NEVER touch or handle the power supply on this instrument due to the high voltage!**

The protection is likely to be impaired if, for example, the apparatus

- ▲Shows visible damage
- ▲Fails to perform the intended measurements
- ▲Has been subjected to prolonged storage under unfavorable conditions
- ▲Has been subjected to severe transport stresses

## **5, Performance**

To ensure that the instrument is working within its specification, especially when making measurements of an important nature, carry out performance checks with particular reference to wavelength and absorbance accuracy. Performance checks are detailed in this manual.

## **6, Radio Interference**

For compliance with the EMC standards referred to in the EC Declaration of Conformity, it is necessary that only shielded cables supplied by us are used when connecting the instrument to computers and accessories.

## CHAPTER 2 INTRODUCTION OF THE INSTRUMENT

### 1. Brief Introduction of the Instrument

This instrument is a single beam, general purpose instrument designed to meet the needs of the Conventional Laboratory, This instrument is ideal for various applications, such as: Chemistry, Biochemistry, Petro-chemistry, Environmental Protection, Food and Beverage Labs, Water and Waste Water Labs and other fields of quality control and research.

This instrument incorporates a 128×64 dots matrix LCD display for photometric results, easy operation and wavelength range of 190nm to 1100nm. This instrument is ideal for measurements in the visible and ultraviolet wavelength region of the electromagnetic spectrum.



Fig 2-1 Front View

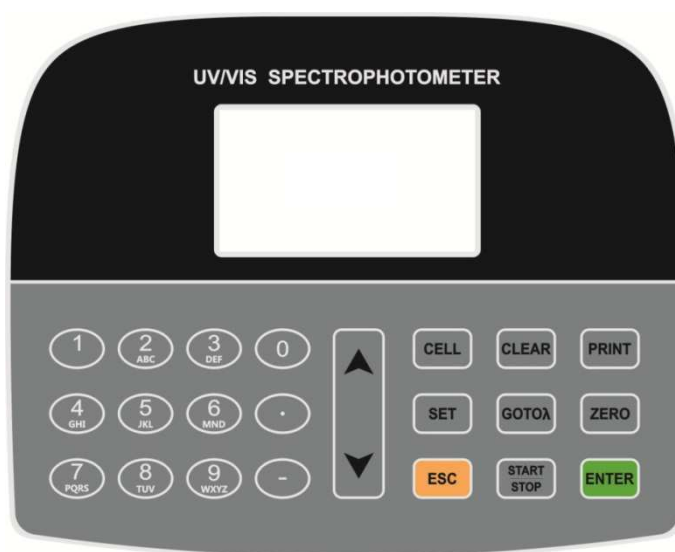


Fig 2-2 Control Panel

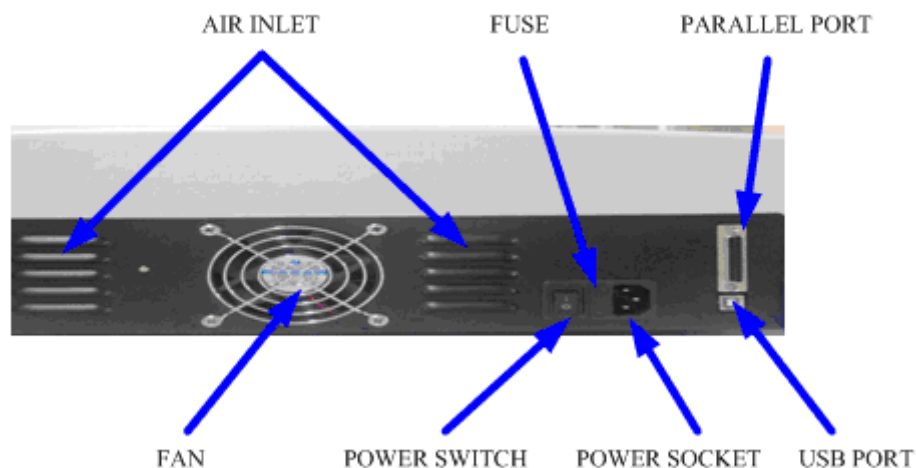


Fig 2-3 Back View

## 2. Working Principle

Different matter has different but special absorbance wavelength point. Also, when at the fixed wavelength point, the absorbance has some relation to the substance's (Always transparent Solution) concentration and its thickness. The relation can be concluded as the following Formula which is called Lambert-Beer Law.

$$T = I/I_0$$

$$A = KCL = -\log I/I_0$$

A Absorbance

C Concentration of the Solution

K Absorbance Coefficient of the Solution

L The length of the Solution in the light path

I The intensity of the light focused on the A/D after it permeate the solution to be measured.

$I_0$  The intensity of the light focused on the A/D after it permeate the Solution.

**Note:** When test, the solvent is usually taken as the Reference Solution and its Transmittance is considered as 100%T. While the Transmittance of the sample to be tested is a relative value which is got comparing to that of the Reference.

## 3. Structure

The spectrophotometer consists of five parts:

- 1) Halogen or deuterium lamps to supply the light;
- 2) A Monochromator to isolate the wavelength of interest and eliminate the unwanted second order radiation;
- 3) A sample compartment to accommodate the sample solution;
- 4) Detector to receive the transmitted light and convert it to an electrical signal; and
- 5) A digital display to indicate absorbance or transmittance. The block diagram (Fig 2.4) below illustrates the relationship between these parts.

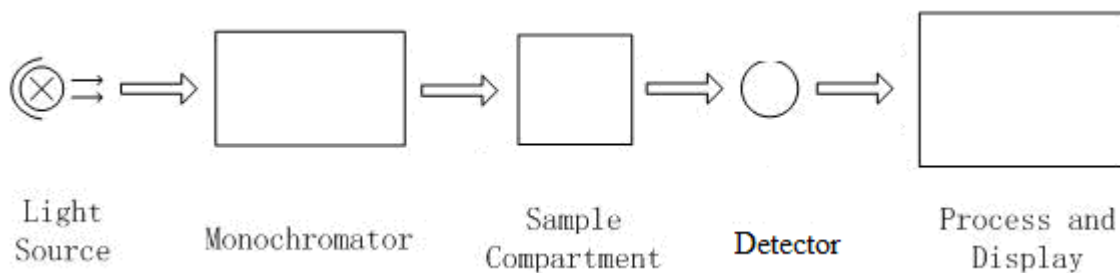


Fig 2.4 Block diagram for the Spectrophotometer

In your spectrophotometer, light from the lamp is focused on the entrance slit of the monochromator where the collimating mirror directs the beam onto the grating. The grating disperses the light beam to produce the spectrum, a portion of which is focused on the exit slit of the monochromator by a collimating mirror. From here the beam is passed to a sample compartment through one of the filters, which helps to eliminate unwanted second order radiation from the diffraction grating. Upon leaving the sample compartment, the beam is passed to the silicon photodiode detector and causes the detector to produce an electrical signal that is displayed on the digital display.

#### 4. Unpacking Instructions

Carefully unpack the contents and check the materials against the following packing list to ensure that you have received everything in good condition.

#### Packing List

Description	Quantity
<i>Spectrophotometer</i> .....	1
<i>Power Cord</i> .....	1
<i>Glass Cuvettes</i> .....	1 Set of 4
<i>Quartz Cuvettes</i> .....	1 Set of 2
<i>Operation Manual</i> .....	1
<i>Software Kit</i> .....	1 Set

*Note: The Printer and Auto-cell Holder mentioned in this manual are all optional accessories, they do not come standard with the instrument.*



## 5. Specifications

Wavelength Range	190-1100nm
Band Width	2nm
Wavelength Accuracy	±0.8nm
Wavelength Repeatability	0.3nm
Wavelength Setting	Auto
Photometric Accuracy	±0.3%T
Photometric Repeatability	0.2%T
Photometric Display Range	0-200%T, -0.3-3.0A
Stability	0.002A/h @ 500nm
Stray Light	≤0.05%T @220nm, 360nm
Data Output Port	USB
Printer Port	Parallel Port
Display	128*64 Dots LCD
Lamps	Deuterium Lamp & Tungsten Halogen Lamp
Detector	Silicon Photodiode
Power Requirement	220V/50Hz or 110V/60Hz
Dimension (L*W*H)	460*360*225mm
Net Weight	18kg
Model UV 4600	Code 6.4200.00

# CHAPTER 3 INSTALL INSTRUMENT

## 1. Environment Required

To ensure the best performance, the following conditions are required:

- The best working temperature range is 16-30°C and the humidity is 45-80%.
- Keep it as far as possible away from the strong magnetic or electrical fields or any electrical device that may generate high-frequency fields.
- Set the unit up in an area that is free of dust, corrosive gases and strong vibrations.
- Remove any obstructions or materials that could hinder the flow of air under and around the instrument.
- The power requirement is 220±22V@50±1Hz or 110±11V@60±1 Hz.
- Use the appropriate power cord and plug into a grounded outlet.
- If the local voltage is not stable enough, a manostat is required.
- Be away from direct sunlight

## 2. Voltage Confirmation



**Be sure to set the instrument' s Voltage Switch at your local power supply, or severe damage may occur! The Voltage Switch is on the bottom of the instrument.**

eg: If your local Voltage is 220V, you should set the Voltage Switch at the following status.(Fig.3-1)



Fig.3-1

If your local Voltage is 110V, you should stir the switch to the Left way by finger, so that it can be set at 110V. (Fig.3-2)



Fig.3-2

### **3. Installation**

#### **Step 1: Check the packing list**

Unpack the contents, check the materials with the Packing List. Any damage or Lost found, please contact us or the local dealer.

#### **Step 2: Position**

Place the instrument on the stable table carefully.

#### **Step 3: Install printer (Optional)**

Make sure the printer is power off; link the printer' s data cable to the instrument' s parallel port.

#### **Step 4: Link the power cord**

Make sure the instrument' s power switch is in the Off condition, link the power cord to the instrument and insert another end in the socket provided with a protective earth contact.

#### **Step 5: Switch on the power**

Check again, make sure that all the links are right. Switch on the power. The instrument can be used to measure after self-test and pre-warm.

# CHAPTER 4 OPERATION INTRODUCTION

## 1. Software Structure

### 1.1 Menu Tree

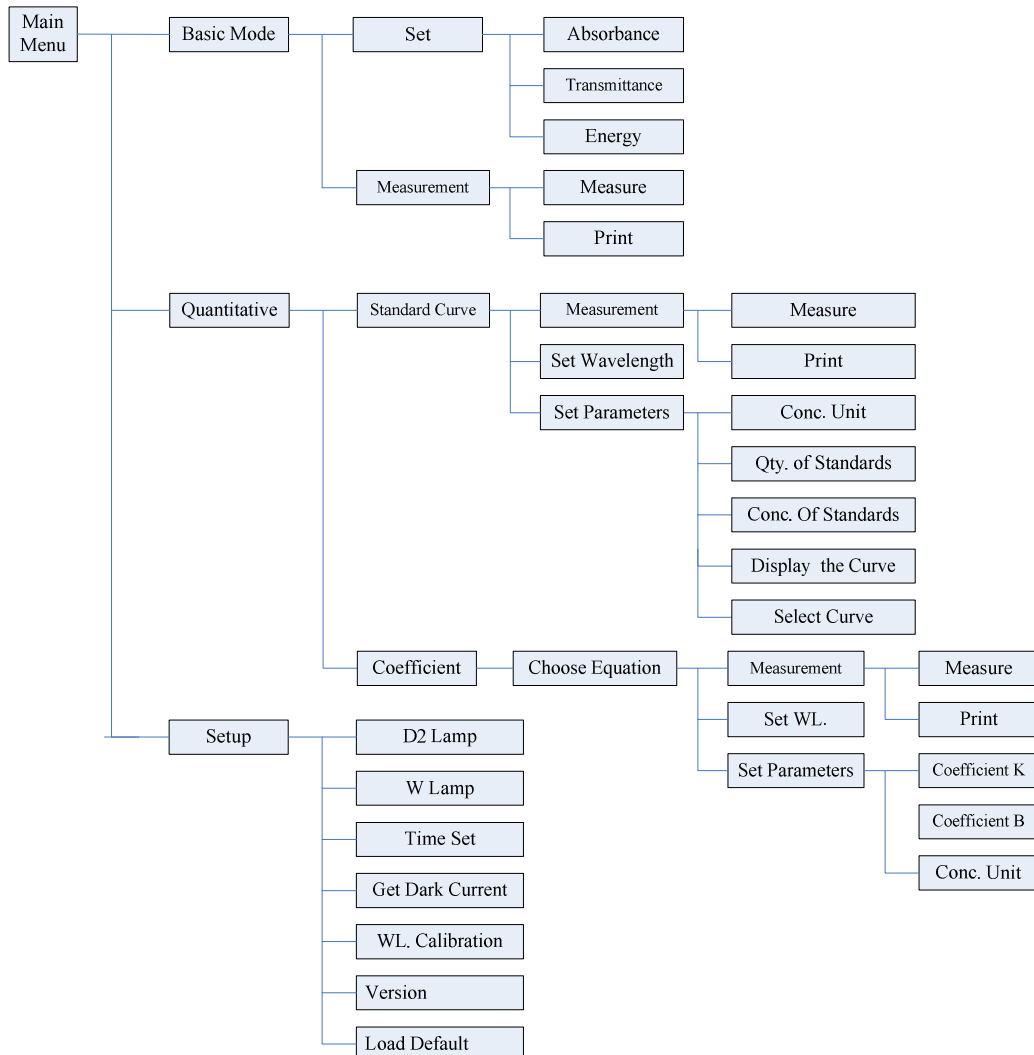


Fig 4-1 Menu Tree

### 1.2 Firmware Functions

The firmware consists of 3 functions: Basic Mode, Quantitative, System Functions.

#### 1) Basic Mode

Measure the Absorbance, Transmittance or Energy in a fixed wavelength point. The test result can be stored in the RAM.

#### 2) Quantitative

**Standard Curve Method:** Set up standard curve by standard samples; use the new curve to measure the unknown sample's concentration. The curves and the test results can be stored in the RAM.


**Coefficient Method:** Input the values of the curve equation' s coefficient, and then measure the unknown sample' s concentration.

### 3) Setup


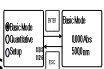
Light source management; Get Dark Current; Calibrate Wavelength and Set time, etc.

## 2. Basic Operation


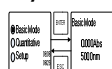
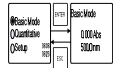
### 2.1 Choose Photometric Mode

Press Arrow key to choose and  to confirm.

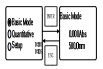
### 2.2 Set Wavelength

In the measurement Interface, press  to set wavelength, use numeric keyboard to input the value, press  to confirm and set 100%T/0Abs automatically.


### 2.3 Set Parameters

Press  to go into the parameter setting interface, press Arrow Key to choose the corresponding menu and use numeric key to input the parameter, press  to confirm, press  to return.

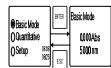
### 2.4 Set the position of auto-cell holder. (The Auto Cell Holder is optional)

In the measurement Interface, press  first, then press the numeric key of 1-8, and then the corresponding cell will be positioned at the light path.

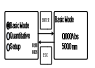
### 2.5 Delete the Input Value

Press  to delete the last character of the input.


### 2.6 Delete the test result and stored data

In the measurement Interface, press  to delete the test result and stored data.

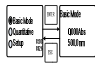
### 2.7 Calibrate 100%T/0Abs

Put the reference in the light path, press  to calibrate 100%T/0Abs.

### 2.8 Measure the samples

Put the samples in the light path and press  to measure.

## 2.9 Print the test result

In the Measurement Interface, press  to print the result..

## 3. Preparation before test

### 3.1 Switch on the instrument and begin the system self-test(Diagnostic Program)

Make sure to remove all the blocks in the light path; Close the compartment lid and switch on the power; Then the system will begin self-test.

**Note: During the Diagnostic Program runs, please DON'T open the lid of the compartment!**

### 3.2 Pre-warm

When the diagnostic program finished, it goes into pre-warm condition. 20 minutes is required before measure.


### 3.3 Check the Cuvettes

The cuvette should be clear and there' s no solution remains on the transparent surface. If the wavelength is below 340nm, **Silicon Cuvette** is a must.

## 4. Measure

### 4.1 Basic Mode

#### Step 1, Go into Basic Mode interface

In the main menu, use Arrow Key to choose "Basic Mode", press  to go into corresponding setting interface. (Fig. 4-2) ;

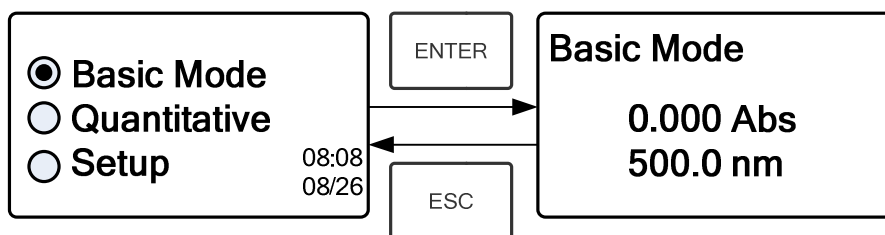
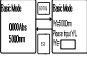



Fig. 4-2

#### Step 2, Set Wavelength

Press  to set Wavelength, input the wavelength value by the numeric Keypad, then press  to confirm. (Fig. 4-3) ;

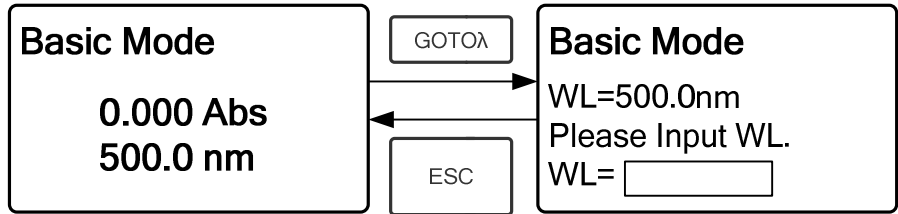


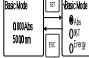


Fig. 4-3

### Step 3, Set Photometric Mode

Press  to go into Parameter Setting Interface, press Array Key to select "Abs", "%T" or "Energy" mode, press  to confirm. Press  to return (Fig. 4-4) ;

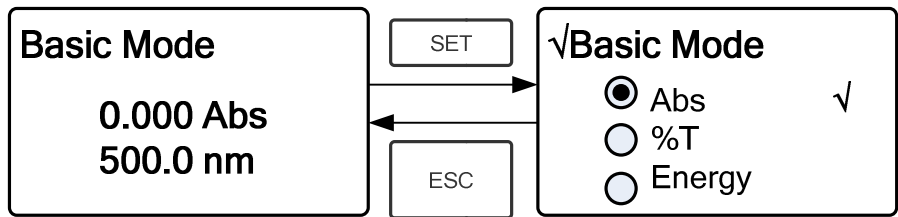



Fig. 4-4

### Step 4, Go into Continuous Measurement Interface

Press  to go into continuous measurement interface (Fig. 4-5) ;

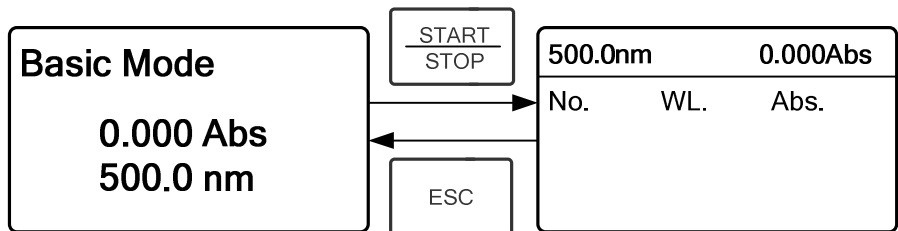


Fig. 4-5

### Step 5, Calibrate 100%T/0Abs

Put the Reference Solution in the light path, press  to calibrate 100%T/0Abs;

## Step 6, Measure Sample

Put the unknown concentration sample in the light path, press  to measure, and the result displays in the data sheet. At the same time, the test result will be stored in the RAM automatically. Repeat this step to finish all the samples measurement.

## Step 7, Print the data

Press  to print the test result.

## Step 8, Delete the data

Press  to delete the test data.

## 4.2 Quantitative

### Step 1: Go into Quantitative Interface

In the main menu, press Arrow Key to choose "Quantitative"; Press  to enter the Quantitative method selecting Interface. (Fig. 4-6) ;

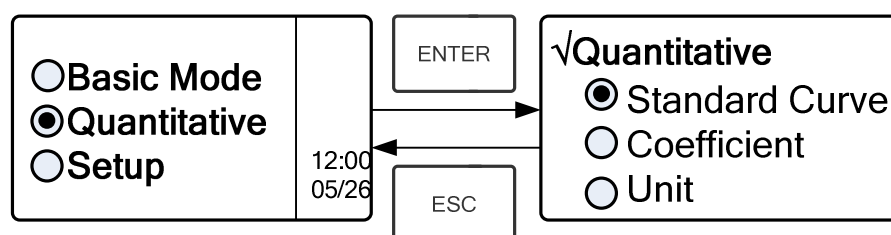


Fig. 4-6

### Step 2: Choose Unit

Move the cursor on "Unit", press  to go into concentration unit choosing interface. (Fig. 4-7) Use Arrow Key to move the cursor on the unit you want, and press

to confirm.



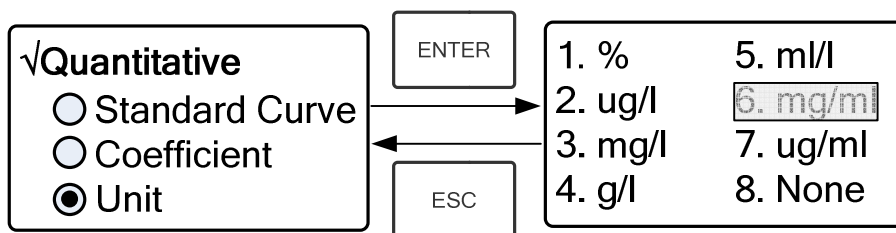



Fig. 4-7

### Step 3: Choose Methods

There are two methods to be selected: Standard Curve and Coefficient.

#### I. Standard Curve:

##### (1) Go into “Standard Curve “Interface

In “Quantitative” menu, press Arrow Key to choose “Standard Curve”; Press  to go into sub menu. (Fig. 4-8) ; Here you can create a new curve, or you can load a saved curve. If you want to delete a saved curve, just choose “Delete Curve” .

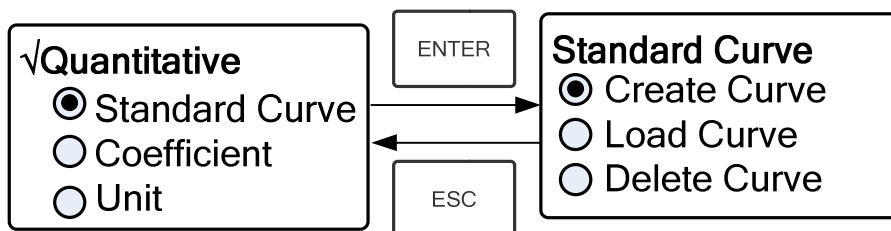



Fig. 4-8

##### (2) Set Wavelength

★ Move the cursor on “Create Curve” , the press  to go into the sample quantity setting interface. (Fig. 4-9)

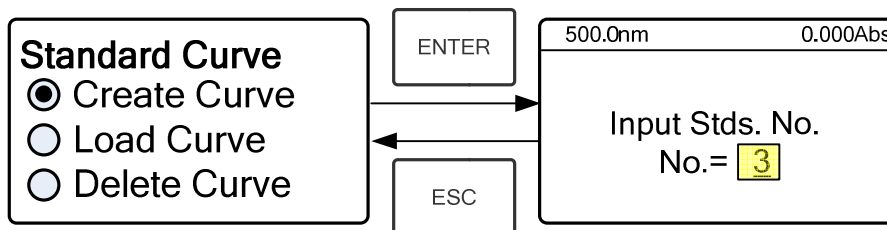




Fig. 4-9

★ Press  to go into the wavelength setting interface, input the WL. Value by the

numeric Keypad and press  to confirm. (Fig. 4-10) ;

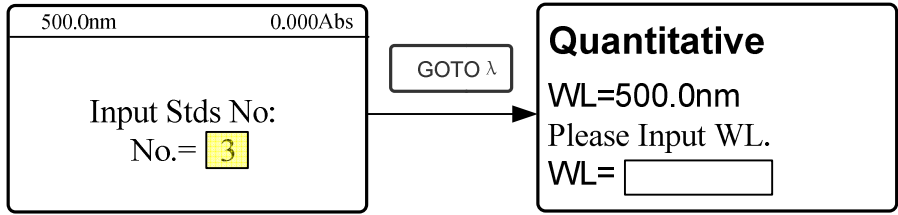
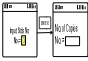


Fig. 4-10

★ Use the numeric keypad or Arrow Key to input the quantity of the standard samples(eg. There are 6 Standard Samples), then press  to confirm. The system will go into the following interface. (Fig. 4-11)

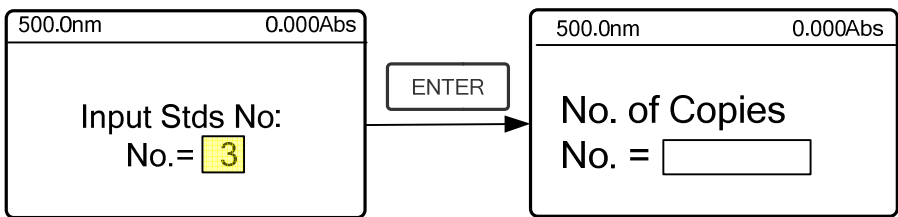



Fig. 4-11

★Use the numeric keypad or Arrow Key to input the test times of each standard sample, then press  to confirm. Then the system will go into the following interface. (Fig. 4-12)

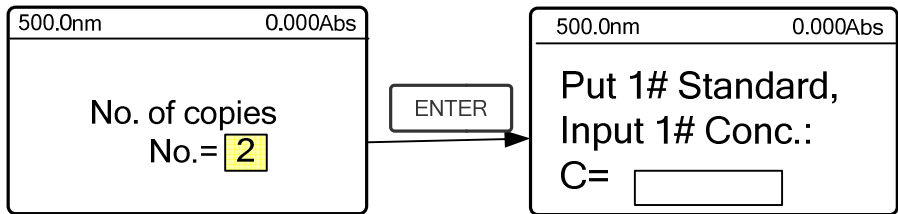




Fig. 4-12

**(3) Calibrate 100%T/0Abs**

Put the Reference Solution in the light path, then press  to calibrate 100%T/0Abs;

**(4) Input Standard Samples' Concentration**

★ Pull the 1# Standard Sample in the light path and input the value of its concentration by the numeric keypad, then press  to confirm. The system will go into the next interface.(Fig. 4-13-1)

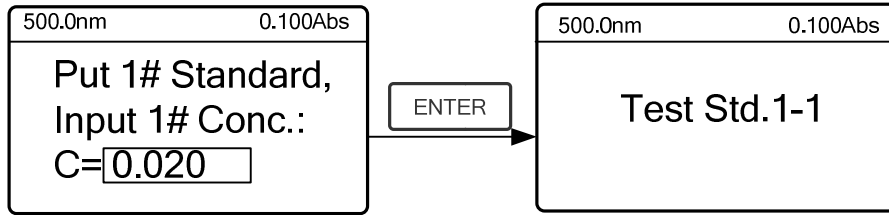


Fig. 4-13-1

★ Press  to measure 1# Standard Sample for the first time, then the system will ask you to test Sample 1# again. (Fig. 4-13-2)

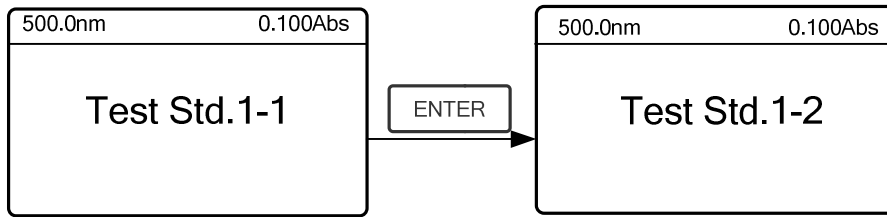


Fig. 4-13-2

★ Press  to measure 1# Standard Sample for the second time. Then it goes into the next Interface. (Fig. 4-14)

*Note: The system will take the Average value of the two results as No.1 Standard Sample' s Absorbance Value.*

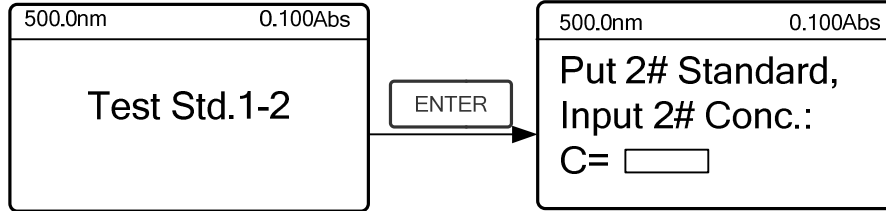
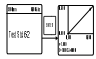
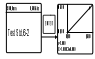


Fig. 4-14

★ Pull the 2# Standard Sample in the light path and input its concentration, then press  to confirm.

★ Measure the other standard samples in the same way. When the last sample is finished, press  to confirm. Then the Standard Curve and its equation will display on the screen automatically. (Fig. 4-15) At the same time, the Curve Equation will be saved in the RAM automatically.

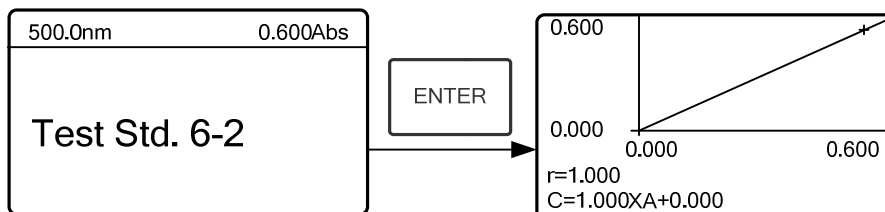



Fig. 4-15

Note: If there's some mistake during the operation, the system will beep 3 times and return the initial interface automatically. Of course, the curve can not be displayed on the screen.

### (5) Measure Samples

★ Pull the Reference solution in the light path and press  to go into the continuous measurement interface. (Fig. 4-16)

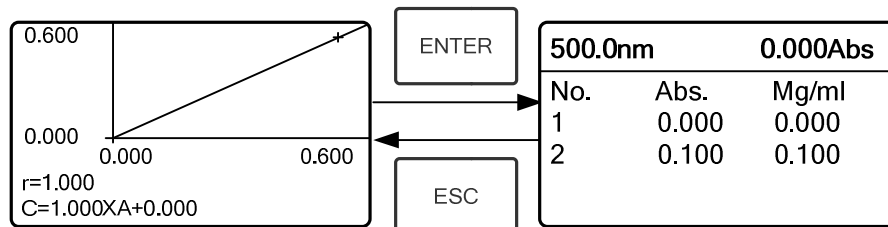




Fig. 4-16

★ Pull the unknown sample in the light path, then press  to measure. The result will be displayed on the screen one by one. At the same time, the data will be saved in the RAM automatically. Total 200 groups of data can be saved.

### (6) Load Curve

All the curve equations will be saved in the RAM automatically. If you want to reload the saved curve, just follow the following steps:

★ Move the cursor on "Load Curve" and press  to confirm. Then the system will go into the curve equation record interface.(Fig. 4-17)

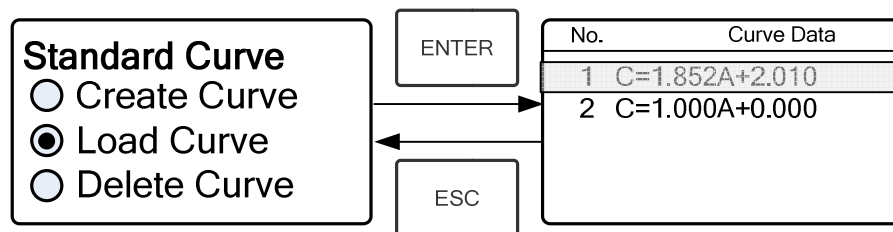



Fig. 4-17

★ Move the cursor on the equation that you need and press  to confirm. Then the corresponding curve will display on the screen. (Fig. 4-18)

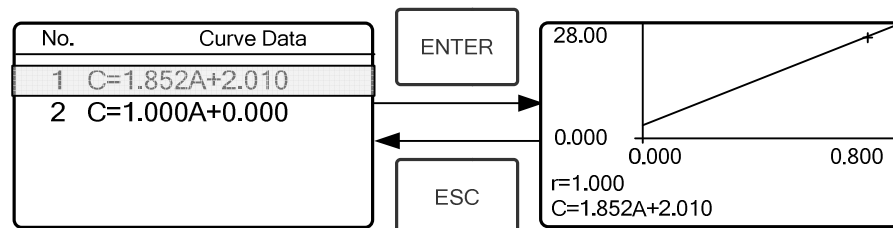



Fig. 4-18

★ Press  to go into sample continuous measure interface and measure samples.

Note: When you reload the curves, the wavelength will be set at the point where you create the Standard Curve automatically.

### (7) Delete Curve

★ Move the cursor on "Delete Curve" and press  to confirm. Then the system goes into the curve equation choosing interface. (Fig. 4-19)

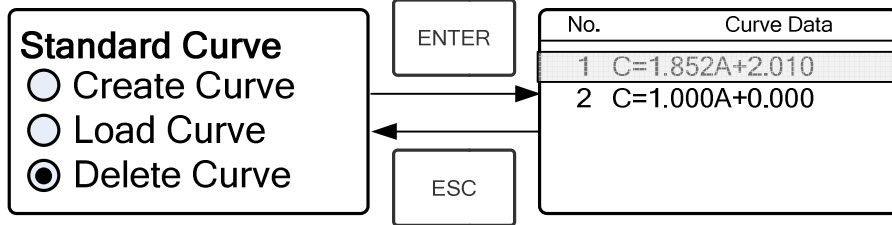



Fig. 4-19

★ Move the cursor on the curve equation that you don't need again and

press , then system will ask you to confirm again.(Fig. 4-20)

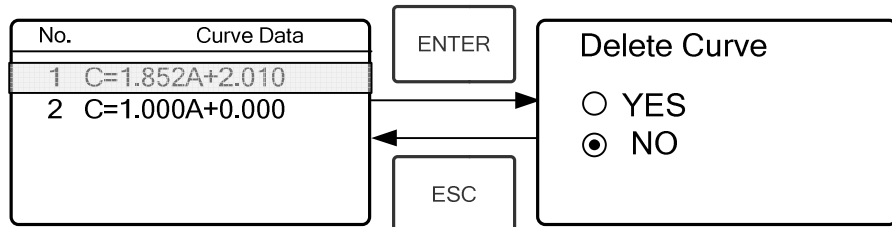






Fig. 4-20

★ Move the cursor on "Yes" and press  to confirm, then the curve equation will be deleted. If you don't want to delete it, just choose "NO" or press  to return.

## II Coefficient Method:

### (1) Go into "Coefficient" method.

Press Arrow Key to choose "Coefficient", press  to go into pre-test interface, then press  to go into the parameter setting interface. (Fig. 4-21) ;

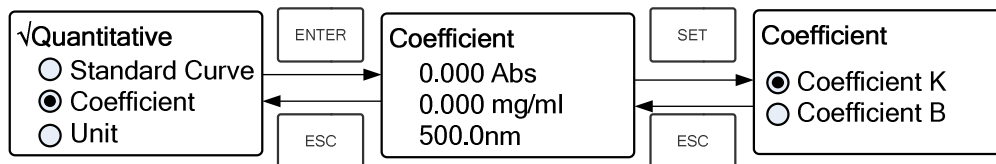




Fig. 4-21

### (2) Parameter Setting

★ Use Arrow Key to move the cursor on "Coefficient K" and press  to go into K setting interface. (Fig. 4-22) ; Input K and press  to confirm, the system will return the former interface automatically.

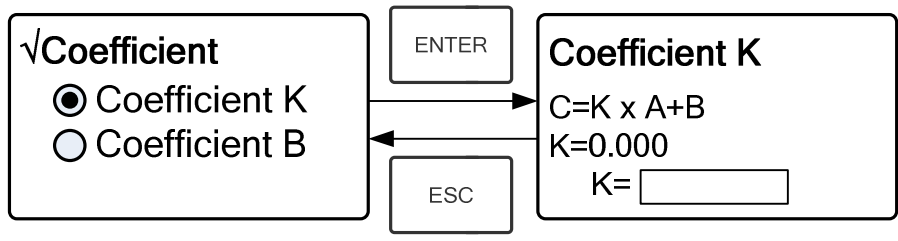





Fig. 4-22

★ Input the value of B in the same way, press  to go back to the pre-test interface.

**(3) Set Wavelength**

Press  to go into wavelength setting interface, input the wavelength value by the numeric keypad and press  to confirm. (Fig. 4-23)

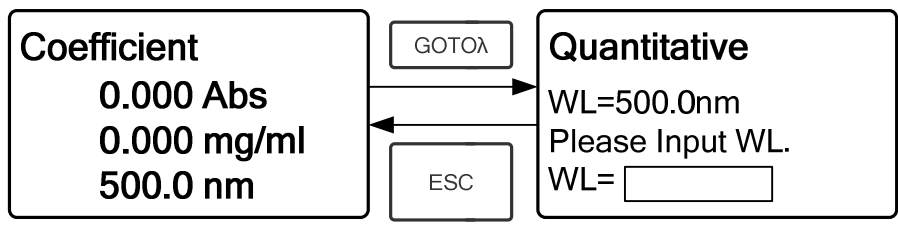

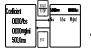


Fig. 4-23

**(4) Calibrate 100%T/0Abs**

Press  to go into continuous measurement interface (Fig. 4-24). Pull the Reference Solution in the light path, then press  to calibrate 100%T/0Abs;

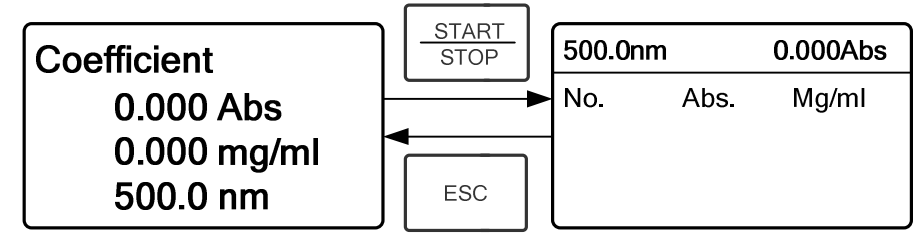



Fig. 4-

24


**(5) Measurement**

Place the samples to be measured in the light path, press  to measure, then the result will display on the screen. Repeat this step to measure other samples. All the test results will be listed on the screen and stored in the RAM automatically. They can be reload even after a sudden power cut.


**Step 4: Print the test result**

In the continuous measurement interface, press  to print the test result.

### Step 5: Delete the test data

In the continuous measurement interface, press  to delete the test result.

## 4.3 SETUP

In the main menu, press Arrow Key to choose "Setup" and press  to go into utility setting interface (Fig. 4-25)

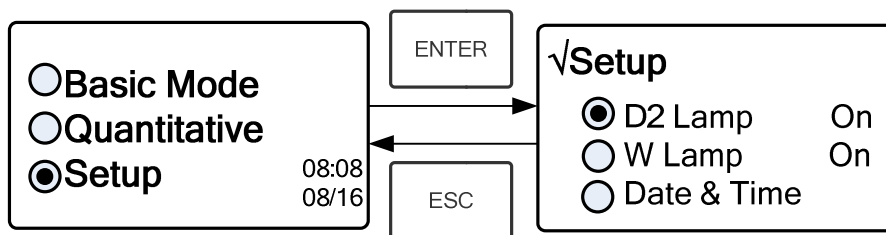

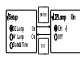
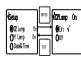


Fig. 4-25

### 1) Power on/off D2 lamp



**When the wavelength point needed is in the range of 340-1100nm, the D2 lamp can be switched off to prolong its life time.**

Use Arrow key to choose "D2 Lamp" and press  to go into D2 lamp setting interface (Fig. 4-26) . Press Arrow key to choose "On " or "Off" and press  to confirm. Press  to return.

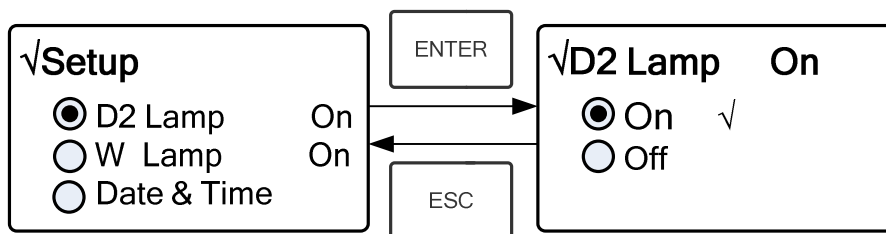





Fig. 4-26

### 2) Power on/off W Lamp



**When the wavelength point needed is in the range of 190-339nm, the W lamp can be switched off to prolong its life time.**

Use the key of Arrow Key to choose "W Lamp", press  to go into W lamp setting interface (Fig. 4-27) . Use the key of Arrow Key to choose "On" or "Off" and press  to confirm. Press  to return.

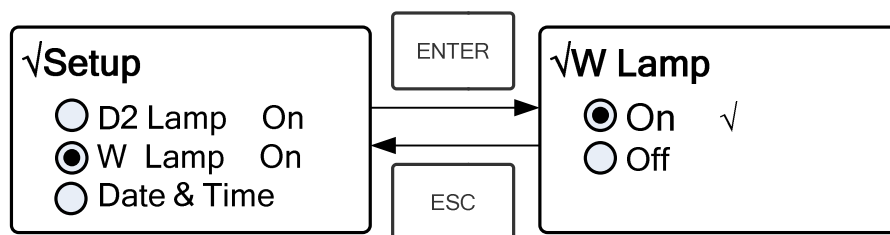




Fig. 4-27

### 3) Set Date & Time

Press Arrow Key to choose "Date & Time" and press  to go into setting interface. (Fig. 4-28) Use Arrow Key to choose setting item and input the value by numeric keypad. (Hour: 0-23, Minute: 0-59, Second: 0-59, Year: 0000-9999, Month: 1-12, Day: 1-31, Week: 1-7) , press  to return.

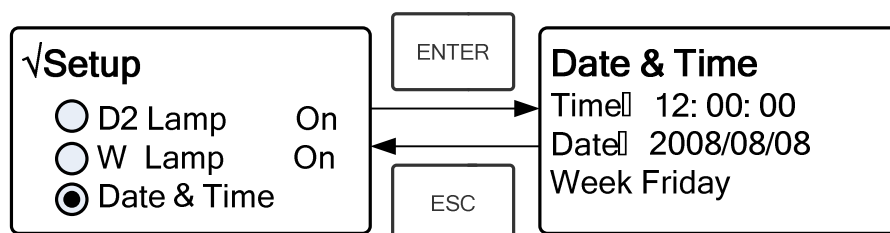



Fig. 4-28

### 4) Calibrate Dark Current



**When the working condition changes, A dark current calibration is required before any measurement.**

Use Arrow key to choose "Dark Current" and press  to confirm. (Fig. 4-29)



**Note: Remove the cuvettes from the light path before this action, don't open the lid of the compartment during the course of calibration.**

Press  to return when it finishes.

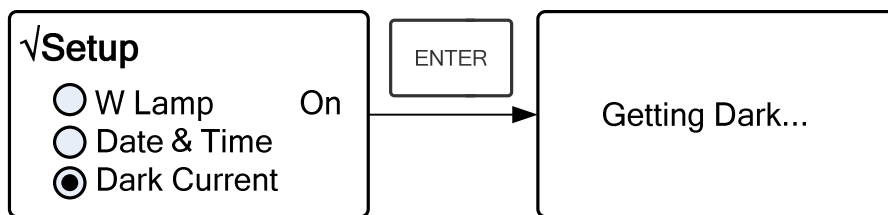



Fig. 4-29

## 5) Wavelength Calibration



**After some time, the energy of lights would drop off and there will be some but slight influence on the test result. In this case, users can reset the wavelength to compensate. We suggest the users to reset the wavelength every one or two months.**

Remove all the blocks in the light path and close the lid of the compartment, press Arrow Key to choose "Calibrate WL." and press  to begin the calibration (Fig. 4-30). The lid of the compartment should not be opened during the calibration. It returns automatically after calibration.

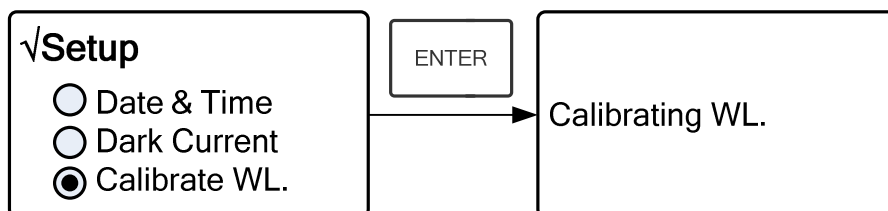




Fig. 4-30



## 6) Lamps Change (Input the lamps changing wavelength point)

**The instrument permits the users to set the lamps changing wavelength point. In the range of 300-400nm, users can choose the lamp freely.**

Press Arrow Key to choose "Lamp Change" and press  to go into corresponding setting interface (Fig.4-31). Input the value of lamps changing wavelength point by the numeric keypad (300—400nm), then press  to confirm and return.

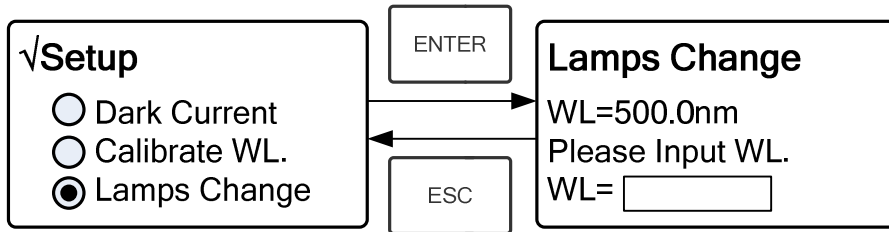




Fig. 4-31

## 7) Load Default

Press Arrow Key to choose "Load Default" and press  to confirm. (Fig. 4-32) ,

press Arrow Key to choose "Yes" and press  to confirm.

Then the system will recover the default setting when leave factory. Choose

"No" with  pressed to cancel the operation and return.

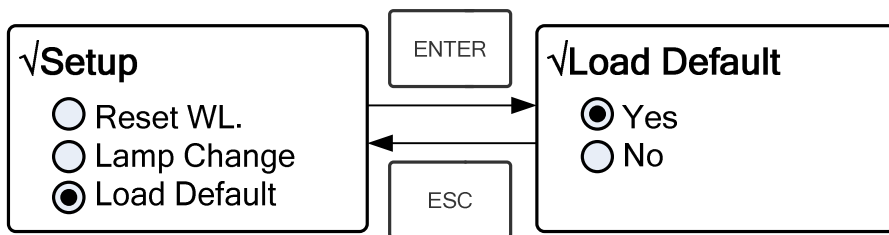




Fig. 4-32

## 8) Version

Press Arrow Key to choose "Version" and press  to confirm. You'll

see the version of the software and hardware. (Fig.4-33). Press  to return.

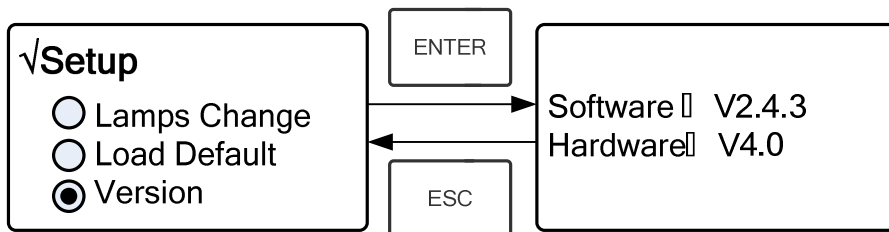


Fig. 4-33

*Note: For the reason of updating, the Version you see on the instrument may be different from the one listed in Fig.4-33.*

# CHAPTER 5 Instrument Maintenance

To keep the instrument work in good condition, constant maintain is needed.

## 1. Daily Maintain

### 1) Check the compartment

After measurement, the cuvettes with sample solutions should be taken out of the compartment in time. Or the volatilization of the solution would make the mirror go moldy. Users must pay more attention to the corrosive sample and liquid easy to volatilize. Any solution remains in the compartment should be wiped off immediately.

### 2) Surface Clean

The cover of the instrument is with paint. Please use wet towel to wipe off the drip on the surface immediately. Organic solution is forbidden to be used to clean the cover. Please wipe off the dirt on the cover timely.

### 3) Clean the cuvettes

After every test or after a solution change, the cuvettes should be cleaned carefully, or the remains on the surface would cause measuring error.

## 2. Troubleshooting

### 1) Dark Current Error when self-test

Possible Cause	Solution
▪ Open the lid of the compartment during the course of self-test.	Close the lid of the compartment and switch on the power again.

### 2) No response after power on

Possible Cause	Solution
▪ Bad contact in power supply	Improve the contact
▪ Fuse melt	Replace a new fuse

### 3) Printer not work, printer error

Possible Cause	Solution
▪ No power supply	Switch on the power supply
▪ Bad contact in power supply	Improve the contact
▪ Bad contact of the data cable	Improve the contact

### 4) No Stable Reading

Possible Cause	Solution
▪ No enough pre-warm	Increase the pre-warm time
▪ Glass cuvettes used in UV Range	Use Silicon Cuvettes.

- |                                        |                             |
|----------------------------------------|-----------------------------|
| ▪ No stable Sample                     | Improve the sample          |
| ▪ Much higher sample concentration     | Dilute the sample           |
| ▪ Low voltage or unstable power supply | Improve the power condition |
| ▪ Lights defect lamp                   | Replace a new               |
| ▪ Light used up                        | Replace a new lamp          |

### 5) Worse Repeatability

#### Possible Reason

- Unstable sample
- Cuvettes polluted

#### Solution

- Improve the sample
- Clean the cuvettes

### 6) Incorrect reading

#### Possible Reason

- Dark Current Error
- Worse matching of the cuvettes

#### Solution

- Re-get the dark current
- Improve the matching of the cuvettes

## 3. Spare part replacement

### Replace the Fuse



**Danger! Be sure to switch off the power and unplug the socket before replacement!**

#### Step 1: Tools preparation

Prepare a 3x75 Flat Blade screwdriver

#### Step2: Switch Off the power supply

Switch off the power supply, and unplug the socket.

#### Step 3: Take out the Fuse Seat

Take out the fuse seat by the screwdriver. (Fig.5-1)



Fig. 5-1

#### Step 4: Replace a new fuse

Pick out the spare fuse and replace it to the working position. (Fig.5-2)

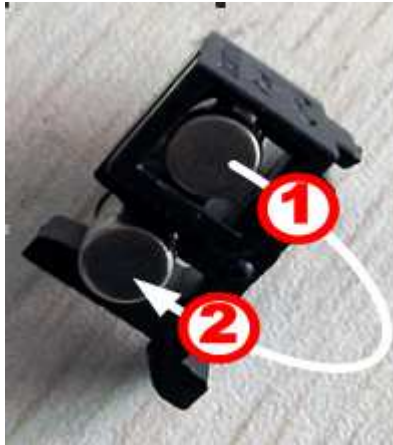


Fig. 5-2

**Step 5: Reset the fuse seat**

Replace the fuse seat in the power socket

**Step 6: Switch on the power**

Plug the socket and switch on the power



Note

Disposal of Waste Electrical and Electronic Equipment by users in private households in the the European Union.

This symbol on the product or on the packing indicates that this cannot be disposed of as household waste. You must dispose of your waste equipment by handing it over to the applicable take-back scheme for the recycling of electrical and electronic equipment. For more information about recycling of this equipment please contact your city office, the shop where you purchased the equipment or your household waste disposal service. The recycling of materials will help to conserve natural resources and ensure that it is recycled in a manner that protects human health and environment.

## Accessories

Code	Articles
6.3200.01	Cuvette holder 1x10mm path light
6.3200.02	Cuvette holder 4x10mm path light
6.3200.03	Cuvette holder 4x50mm path light
6.3200.04	Cuvette holder 4x100mm path light
6.6900.01	Cuvette holder 8x10mm path light, automatic, only for Model UV-6900
6.3200.04	Cuvette holder 1x10mm path light for recirculating water
6.3200.05	Micro Cuvette holder , adjustable
6.3200.06	Transparent Film holder
6.3200.07	Tube holder , max. 20 mm Ø
6.3200.08	Peltier System 15°C to 65°C
6.3200.09	Sipper system 1 to 256s
6.3200.10	Peltier / Sipper system
6.6900.07*	Micro printer
6.3000.01	PC Software-Quantitative/Scanning/Kinetics/Multi- wavelength/AND/Protein
8.9749.00*	PC with monitor TFT 19"
6.6900.03	Halogen lamp visible, all models
6.6900.08	Deuterium lamp, ultraviolet, all models except UV-4000

Code	Articles
6.3200.11	Cuvette with lid, glass, 1.7ml, path light 5mm, 4pcs
6.6900.05	Cuvette with lid, glass, 3.5ml, path light 10mm, 4pcs
6.3200.12	Cuvette with lid, glass, 7ml, path light 20mm, 4pcs
6.3200.13	Cuvette with lid, glass, 10.5ml, path light 30mm, 4pcs
6.3200.15	Cuvette with lid, glass, 17.5ml, path light 50mm, 4pcs
6.3200.16	Cuvette with lid, glass, 34ml, path light 100mm, 4pcs
1.9363.00	Cuvette with lid, plastic, 3.5ml, path light 10mm, 100pcs
6.4000.01	Cuvette with lid, quartz ,1.7ml path light 5mm, 2pcs
6.6900.06	Cuvette with lid, quartz, 3.5ml path light 10mm, 2pcs
6.4000.02	Cuvette with lid, quartz, 7ml path light 20mm, 2pcs
6.4000.03	Cuvette with lid, quartz 10.5ml path light 30mm, 2pcs
6.4000.05	Cuvette with lid, quartz, 17.5ml path light 50mm, 2pcs
6.4000.06	Cuvette with lid, quartz, 34ml path light 100mm, 2pcs
6.4000.07	Cuvette micro 12,5*12,5*45, 100µl, 2pcs
6.4000.08	Cuvette micro 12,5*12,5*45, 200µl, 2pcs
6.4000.10	Cuvette micro 12,5*12,5*45, 700µl, 2pcs

## "CE" DECLARATION OF CONFORMITY

DINTER S. A. *DINKO Instruments* c/ Encarnació, 123-125 / 08024 – Barcelona

Declare that the products mentioned in attached list, to which this declaration relates, comply with the essential safety requirements of the relevant European Directive.

- Low Voltage Directive D2006/95/CEE, of December 12, 2006

- Electromagnetic compatibility (EC), relative to the Directive for electromagnetic compatibility 2004/108/CEE, of December 15, 2004

- Safety for electrical devices of measurement, control and laboratory use. Requirements concerning to CEM. EN 61326

- Safety rules for the electric devices of measurement, control and laboratory use. Part I. General requirements EN 61010-1

Mounting and connecting instructions defined in catalogues and technical data sheets must be respected by the user.

Name: Joan A. Bravo  
Position: Technical Manager

Josep X. Sensada  
Quality Manager

Signature:



Model: Spectrophotometer UV-VIS 4600

## **OTHER DINKO APPARATUS**

- Colorimeters
- Extractor for mince analysis
- Heater Metallic Blocks
- Heater Plates
- Infrared Ovens
- Kits for Water Analysis
- Magnetic Stirrers.
- Muffle Furnaces to 1200°C
- Nephelometers
- Orbital Shakers
- Peristaltic Pumps
- Photometers
- Rotary Stirres
- Sand Baths
- Spectrophotometers
- Stirrers
- Temperature Controllers
- Timers
- Tube Vibrators
- Turbidimeters
- Turn Dishes
- Ultrasonic Baths
- Vacuum Pumps